



CB500 and B600 - How to Change Your Account Password

N. Duay, D. Olson; 28 Jan 2021


- Click the **“Change User”** button in the lower right of **“Icon NMR: Automation”**
- Logon with your directory name and your temporary password: **nmr**
- Once logged on, click **“Options”** at the top of the screen

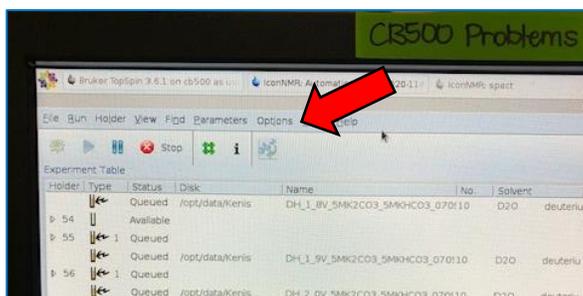

- Then click **“Change Additional User’s Password”**

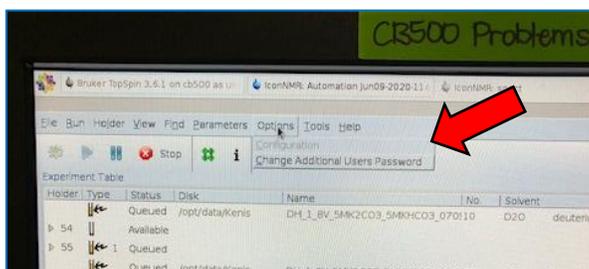

- Enter the current password (**nmr**) and then your new password twice.
- Hit **“Enter”**.

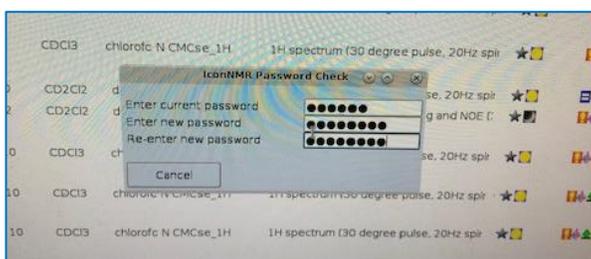

- To check your new password, log out by clicking the **“Change User”** icon then log on as usual with your new password.